


DIGITAL MICROSCOPE INSTALLATION MANUAL

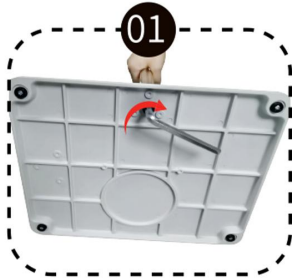

01
Mounting the Column

First loosen the lower nut, then insert the column through the base plate, secure the top and base plate with hex bolts.



02
Installing the Locking Ring

Loosen the locking ring screw, slide to desired height on the column and tighten.



03
Attaching the Focus Wheel

Mount the focus wheel bracket to the designated position on the column.



04
Camera Installation

Remove the protective cap from the camera port. Thread and tighten the lens assembly onto the camera.



05
Locking Mechanism

Loosen the focus bracket screw, adjust position then re-tighten.




06
Magnification Adjustment

Turn left gear for fine magnification, right for reduction.




07
LED Light Assembly

Remove the light cover retention nut. Insert the LED into the slot below the lens and secure.

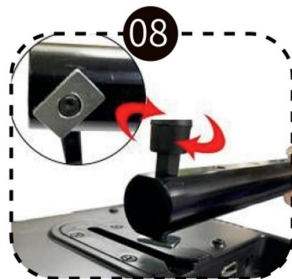

08
Display Unit

Align with the display backplate slot and press into position. Tighten the front panel fastening nut.

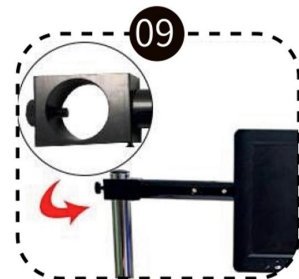

09
Crossbar Mounting

Slide the display-mounted crossbar into the column top groove.

Actual product accessories may vary slightly in appearance from illustrations in this manual.